

Agilent 7000E Triple Quadrupole GC/MS System



Compatible with both Agilent 8890 and Intuvo 9000 GC systems, the Agilent 7000E mass spectrometer (MS) is the best choice for routine GC/MS/MS analysis with a low 4 fg octafluoronaphthalene instrument detection level (IDL) specification. The 4 fg IDL is demonstrated at installation, verifying total system performance of the automatic liquid sampler (ALS), GC, and MS.

Improved mass spec intelligence combined with over five decades of innovation and quality embody the Agilent 7000E GC/TQ as it provides robust, day-in, day-out performance. The 7000E GC/TQ is powered by new technologies that drive maximum lab productivity, so you can focus on moving your lab forward.

7000E triple quadrupole GC/MS

Parameter	Value
Mode of Operation	EI standard, CI optional
Ion Source Material	Noncoated, proprietary inert source
Ion Source Temperature	150 to 350 °C
Filaments	Dual filaments for EI
Source Cleaning	Automated and vent-free with patented Agilent JetClean option
Electron Energy	10 to 300 eV
Mass Filters (2)	Proprietary monolithic hyperbolic gold-coated quadrupole
Mass Axis Stability	<±0.10 u over 24 hours (10 to 40 °C)
Quadrupole Temperature	106 to 200 °C
Mass Range	<i>m/z</i> 10 to 1,050
Dynamic Range	>6.0 ×10 ⁷ counts, resulting in up to seven orders of linear dynamic range from the LOD
Resolution	Selectable, 0.7 to 2.5 Daltons, default tune Settable, 0.4 to 4.0 Daltons, custom tune
Scan Rate	Up to 20,000 u/s
Tuning	Autotune or manual
Detector	Agilent triple-axis HED-EM with extended-life EM and dynamically ramped iris
MRM Speed	800 transitions/sec
Minimum MRM Dwell	0.5 msec
Collision Cell	Linear hexapole
Collision Cell Gas	Nitrogen with helium quench gas
Collision Energy	Selectable up to 60 eV
Vacuum System	Dual stage turbomolecular pump Total gas flow up to 8 mL/min
Software	Agilent MassHunter Acquisition, data handling (quantitative/qualitative) and reporting Pesticides and Environmental Pollutants MRM database with over 8,000 optimized transitions (optional)
Data Acquisition Modes	Full scan, SIM, MRM, dMRM, tMRM, dMRM/scan, and tMRM/scan

8890 GC

For more specifications, refer to the GC data sheet

Parameter	Value
Injector	Split/splitless, multimode inlet, PTV, and others
Autosampler	Agilent 7693 ALS, CTC PAL3, Agilent 7697A headspace sampler
Oven Temperature	Ambient +4 to 450 °C
Oven Ramps/Plateaus	Supports 20 oven ramps and 21 plateaus Negative ramps are allowed
Ramp Rate	120 °C/min (200 +V), 75 °C/min (120 V)
Capillary Flow Technology	Effluent splitting, backflushing, column switching
Retention Time Locking (RTL)	RTL-ready, constant flow or pressure

Intuvo 9000 GC

For more specifications, refer to the GC data sheet

Parameter	Value
Injector	Split/splitless and multimode inlet
Autosampler	Agilent 7693 ALS, CTC PAL3, Agilent 7697A headspace sampler
Oven Temperature	Ambient +10 to 450 °C
Oven Ramps/Plateaus	Supports 20 oven ramps and 21 plateaus Negative ramps are allowed
Ramp Rate	250 °C/min
Capillary Flow Technology	Effluent splitting, backflushing, column switching
Retention Time Locking (RTL)	RTL-ready, constant flow or pressure

Integrated GC/MS system features

Parameter	Value
Early Maintenance Feedback (EMF)	Monitors GC and MS resources: injection counter, operation times, and electronic logs to aid planned maintenance
Parts Finder	Easy, convenient access to pertinent consumables part numbers
Quick Vent	Automated, rapid venting of the MS
Eco-Friendly Operation	User-scheduled sleep/wake mode saves carrier gas and power
Integrated Calculators	Vapor volume calculator, solvent vent calculator, method translator, and so on

Instrument dimensions

Parameter	Value
7000E MS	35 cm (w) × 86 cm (d) × 47 cm (h), Weight: 59 kg Additional space for the data system and optional printer
Mechanical Pump	18 cm (w) × 35 cm (d) × 28 cm (h), Weight: 21.5 kg
8890 GC	58 cm (w) × 51 cm (d) × 49 cm (h), Weight: 49 kg

Reference specifications*

Parameter	Value
EI MRM S/N (He Carrier)	1 µL of 10 fg/µL OFN produces >1,500:1 RMS S/N for the transition of m/z 272 → 222 1 µL of 100 fg/µL of OFN produces >15,000:1 RMS S/N for the transition of m/z 272 → 222
PCI MRM S/N (He Carrier)	1 µL of 100 fg/µL BZP produces >50:1 RMS S/N for the transition of m/z 183 → 105 (CH_4)
NCI SIM S/N (He Carrier)	1 µL of 100 fg/µL OFN produces >2,000:1 RMS S/N for m/z 272 (CH_4)

*Reference specifications represent typical performance, and are not confirmed at installation.

Installation checkout specifications

IDL is a more accurate indication of true sensitivity (minimum detectable quantity) than signal-to-noise (S/N), particularly when background noise levels are very low, as with MS/MS measurements when only standard is injected.

IDL verification is a more extensive (eight consecutive injections versus one, in the case of S/N) and reliable test that is performed at installation to ensure proper system qualification.

Parameter	Value
EI MRM IDL	4 fg or less octafluoronaphthalene (OFN) Statistically derived at 99% confidence level from the area precision of eight sequential splitless injections* of 1 µL, 10 fg/µL OFN standard. MS/MS transition of m/z 272 → 222 100 msec dwell time
HydroInert EI MRM IDL (H_2 carrier)	8 fg or less octafluoronaphthalene (OFN) Statistically derived at 99% confidence level from the area precision of eight sequential splitless injections* of 1 µL, 10 fg/µL OFN standard. MS/MS transition of m/z 272 → 222 100 msec dwell time
PCI MRM S/N	1 µL of 5 pg/µL benzophenone (BZP) produces >2,500:1 RMS S/N for the MS/MS transition of m/z 183 → 105 (CH_4)

–EI MRM IDL and PCI MRM S/N tests were carried out on a 30 m × 0.25 mm, 0.25 µm column.

–HydroInert EI MRM IDL test carried out on a 20 m × 0.18 mm, 0.18 µm column.

*IDL specification is only demonstrated if an autosampler is part of the installed system. If an autosampler is not present, the EI MRM S/N specification is performed.

www.agilent.com/chem/gc-ms-ms

This information is subject to change without notice.

DE62937435

© Agilent Technologies, Inc. 2022
Printed in the USA, June 24, 2022
5994-0481EN